

Peroxidase Block

Cat# NB-23-00192-1 (100 ml)

Cat# NB-23-00192-2 (18 ml)

Intended Use

Peroxidase Block Kit is used to block endogenous peroxidase enzymatic activity which can be used in paraffin-embedded tissue sections. **Peroxidase Block Kit** is provided with the ready-to-use solution in a dropper bottle.

Kit Components

Product	Catalog No.	Size	Number of Slide Tests
Peroxidase Block Kit (RTU)	NB-23-00192-2	18mL	180 slides
Peroxidase Block Kit (RTU)	NB-23-00192-1	100mL	1000 slides

Recommended Protocol

The Peroxidase blocking reagent should be used prior to application of other pretreatments or antibody to tissue.

After deparaffinization:

- 1) Apply 2 drops (50 μ L/ drop) of **Peroxidase Block** to cover the tissue and incubate for 10 minutes at room temperature (25°C). Optimizing block time for your experiments is recommended.
- 2) Rinse well with distilled water.
- 3) Proceed to next step.

Storage

Store at 2-8°C. Return reagent back to 2-8°C after each use.

Precautious

Please wear gloves and take other necessary precautions.

For reference only

For Research Use Only. Not for Diagnostic or Therapeutic Use.